

HIV Ag/Ab Combo ELISA KIT


—Research Use Only—



- HIV 抗原抗体同時測定ができます
- WHOが推奨する第4世代の製品です
- 感染初期の検出が可能です

定価：@55,000円（税別）

HIV-1 p24 Antigen				HIV-1 patient sample			
pg/mL	well 1	well 2	average	希釈	well 1	well 2	average
0	0.046	0.043	0.02967	0	0.046	0.043	0.02967
2	0.069	0.067	0.04533	0.00002	0.131	0.130	0.08700
5	0.150	0.138	0.09600	0.00005	0.257	0.255	0.17067
10	0.189	0.179	0.12267	0.0001	0.474	0.466	0.31333
20	0.339	0.314	0.21767	0.0002	0.839	0.832	0.55700
50	0.796	0.732	0.50933	0.0005	1.736	1.734	1.15667
100	1.461	1.373	0.94467	0.001	2.595	2.557	1.71733
200	2.693	2.557	1.75000	0.002	3.403	3.346	2.24967

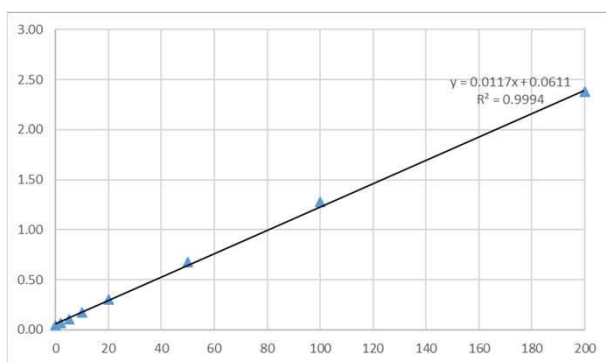

Antigen detection range:
2.0pg/ml～200pg/ml

Rimco

リムコ株式会社

〒904-2234

沖縄県うるま市州崎12-75-106

Tel 098-923-2071 Fax098-923-2072

Email: info@rimco.jp


HIV Ag/Ab Combo ELISA KIT

—Research Use Only—

Results for the Rimco HIV combo ELISA assay in HIV patient and healthy volunteer plasma


A社	Rimco HIV combo	
	Positive results	Negative results
Positive results	34	0
Negative results	0	19
Sensitivity	1.00	
Specificity	1.00	

RIMCO HIV combo vs A社

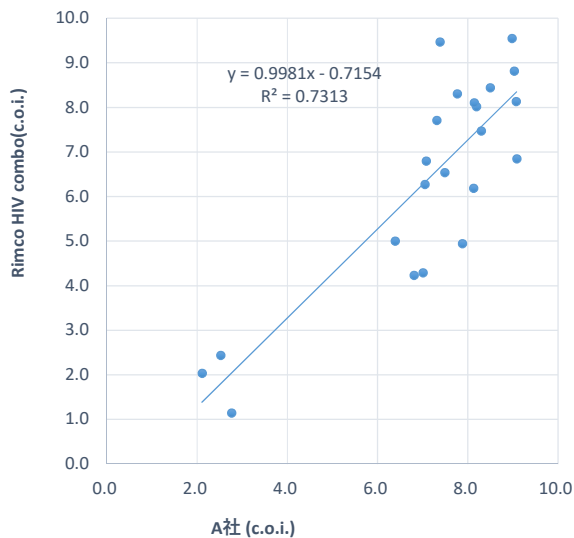

Correlation between the results of the Rimco HIV combo ELISA and A社 for 22 field-collected HIV plasma samples. The correlation ratio(R2) between the Rimco HIV combo ELISA and A社 results was 0.731.

RIMCO HIV combo vs. A社

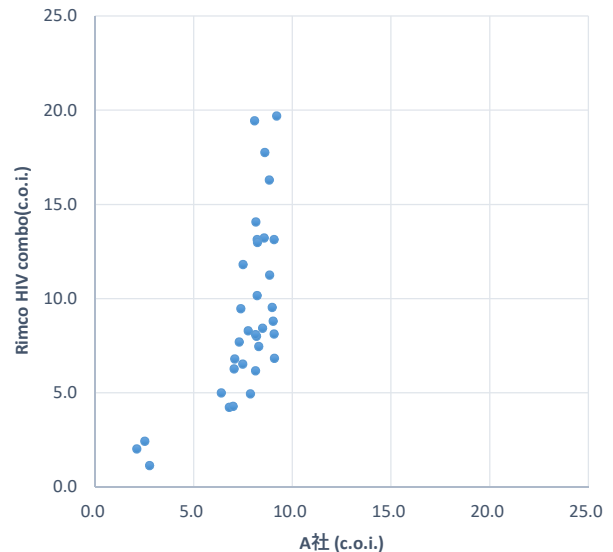

Correlation between the results of the Rimco HIV combo ELISA and A社 for 34 field-collected HIV plasma samples.

Rimco

リムコ株式会社

〒904-2234

沖縄県うるま市州崎12-75-106

Tel 098-923-2071 Fax098-923-2072

Email: info@rimco.jp